

## Cyanoacrylate Gel Preclinical Component Specification

- Convenient, ready-to-use adhesive gel (alternative to cranioplastic/dental cement)
- Adheres Cannula pedestals to the skull of laboratory animals without mounting screws
- Ideal for mice and other small animals when using Short Pedestal Cannulas
- A 20 gram tube supplies enough gel for approximately 100 animals

NOTE: Cannot be used in the same way as cranioplastic/dental cement

### *Advantages:*

- No messy mixing required
- Faster drying time compared to traditional dental cement
- Less weight on the animal's skull minimizes stress and trauma
- Reduced waste
- Economical - more uses per tube

### *Disadvantages:*

- Should not be used with a swivel or other torque inducing equipment
- Cannot be used to build up under the cannula to adjust depth
- If too much is used, curing time is affected
- Shorter adhesive duration than dental cement and screws

PART NUMBER	TYPE	TUBE SIZE
300CAGS	GEL	3 GRAM
300CAG	GEL	20 GRAM